

Standard Cheese Whey Samples – Filter Guide

START



Material for whey filtration:

- Container for taking whey sample
- Vial for filtered sample
- Syringe to extract whey for filtering (10mL)
- Filter 0,45 μm to filter whey sample
- Pen for labeling your sample

1



Disinfect your hands

Fill the syringe with the sample to be filtered

2



Screw the filled syringe on the filter


3



Slowly press down on the syringe to start filtration

5-10mL are needed

SEND

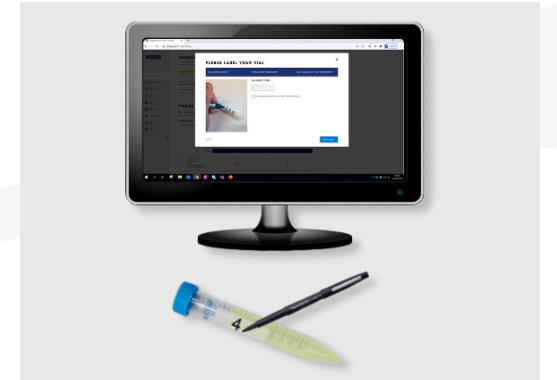

Label the vial with the number you input into PhageWatch™ and send